

Storing and Hiding Data Runs in Logger Pro

- Please do not remove this sheet from the table.

1. Storing data runs in Logger Pro

You can store multiple runs of data collected by Logger Pro in memory for comparison. Normally after each data acquisition run the data are cleared before the next run begins. If you want to keep the data in memory:

- On the main menu in *Logger Pro* click on **Experiment** and then choose **Store Latest Run**.
- The **Store Latest Run** option saves the most recently collected data in memory allowing you to take another data set without losing the first one. It is then possible to compare experiments.
- Runs are stored sequentially as Run 1, Run 2, and so forth.

2. Hiding data runs in Logger Pro

You can hide a stored data run and prevent it from being displayed by *Logger Pro* by clicking on **Hide Data Set** in the **Data** menu and then choosing the specific run from the list that pops open.

- To redisplay a hidden run click on **Show Data Set** in the **Data** menu and choose the run from the list that pops open.

Note: To delete a stored run click on **Delete Data Set** in the **Data** menu and choose the run from the list that pops open.